Transmission Reliability of Implantable Wireless Temperature Loggers for Laboratory Rats

Ásgeir Bjarnason^{1,2}, Hákon Guðmundsson², Sigurður Gunnlaugsson², Andrés Gunnarsson², Kjaran Sveinsson², Sigmar Guðbjörnsson²

1 Tampere University of Technology, Department of Electronics and Communication Engineering, Korkeakoulunkatu 3, Tampere, Finland E-mail: asgeir.bjarnason at tut.fi, sigmar at star-oddi.com 2 Star-Oddi Ltd. Garðabær, Iceland

Objectives


- To measure the transmission reliability of low-power implantable wireless temperature loggers in laboratory rats.
- To test and verify a simple method that enhances the reliability of the wireless transmission.
- To measure the transmission power and pattern of the system.

Methods

Each cage has an antenna placed below it and an RF Every receiver handle up to 10 subjects per cage. The RF box sends information to the PAN.



connected tcomputer via USB. It receives a signal from the RF box attached to each cage. Every PAN can handle up to 64 RF boxes.

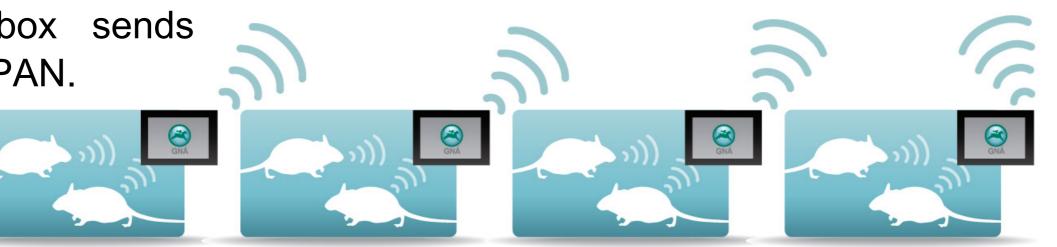

The RF transmission is done through a custom-made protocol that uses a 500kHz output frequency to send the ID, temperature bandwidth and losses due to the data and CRC. In transmission mode the average current consumption is 2.5mA. The receiving antenna placed under the cage in the experiments measured 250x350mm.

The 500kHz frequency band is chosen biological tissue. measurement (n).

since it offers a trade-off between power, To improve the throughput of transmission a simple method was developed, where the three last measurements (n, n-1, n-2) are sent with each transmission instead of a single


Results


Transmission throughput was initially measured at 94% on average ranging from a minimum of 87% to a maximum of 97% transmitting a single measurement (n).

By sending three measurements (n, n-1, n-2) it was possible to increase the reliability of the transmission up to 98.4% on average (min=97.7%, max=98.8%).

The average transmission power inside of the rat's cage was measured with an unamplfied receiver in free space conditions at a -38.53dBm where transmission strongest was inside the receiving antenna's loop -35.5dBm on average.


implanted the Once absolute transmission power decreases slightly. On the other hand the transmission pattern becomes more uniform and is not as vulnerable to orientation of the antenna free in space as conditions.

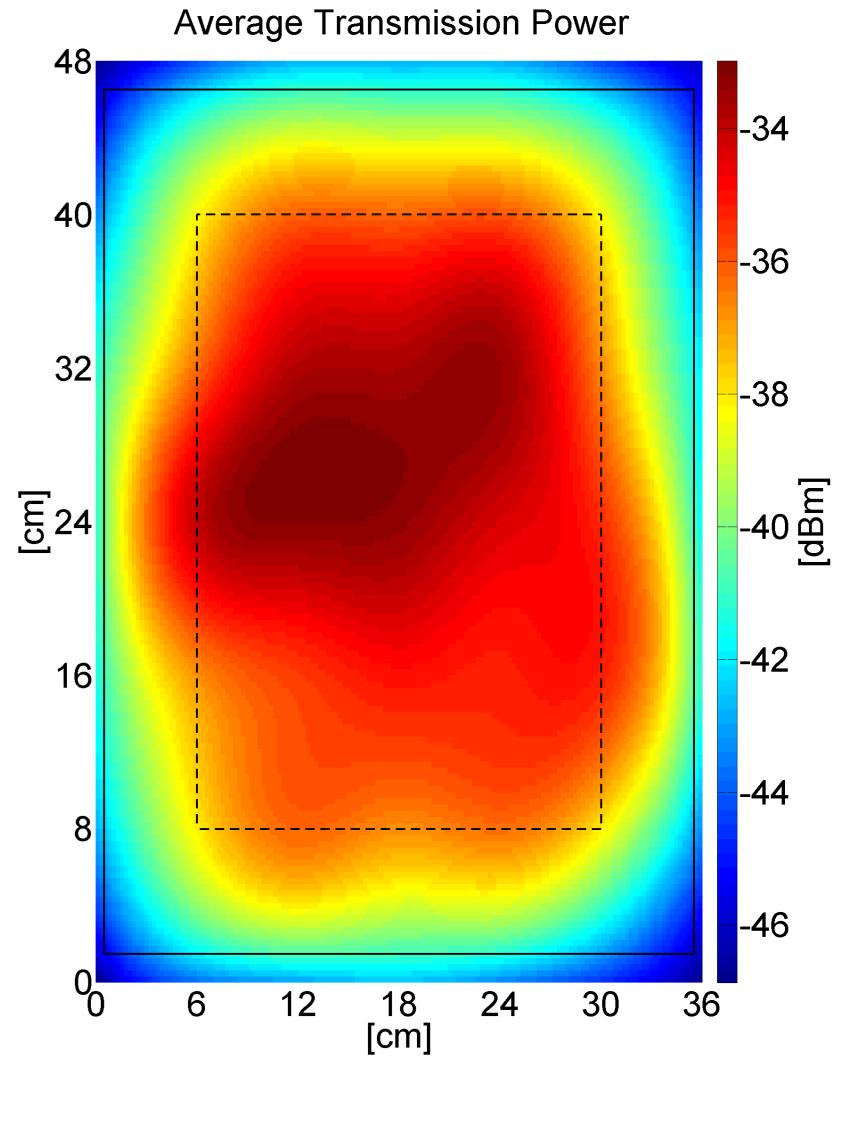

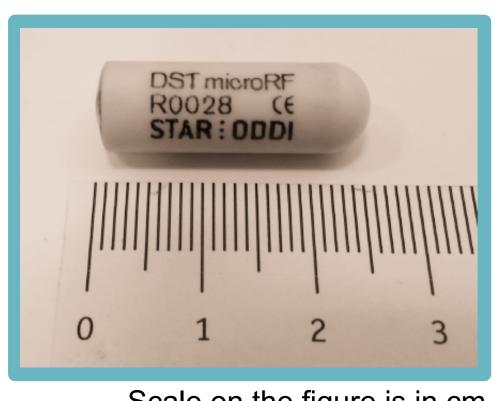



Scale on the figure is in cm

storage tag (DST) is cylindrically a shaped ceramic capsule 8.3x25.4mm) (dimensions: with bio-compatible filled 3.3gr weighs epoxy. including nonrechargeable battery.

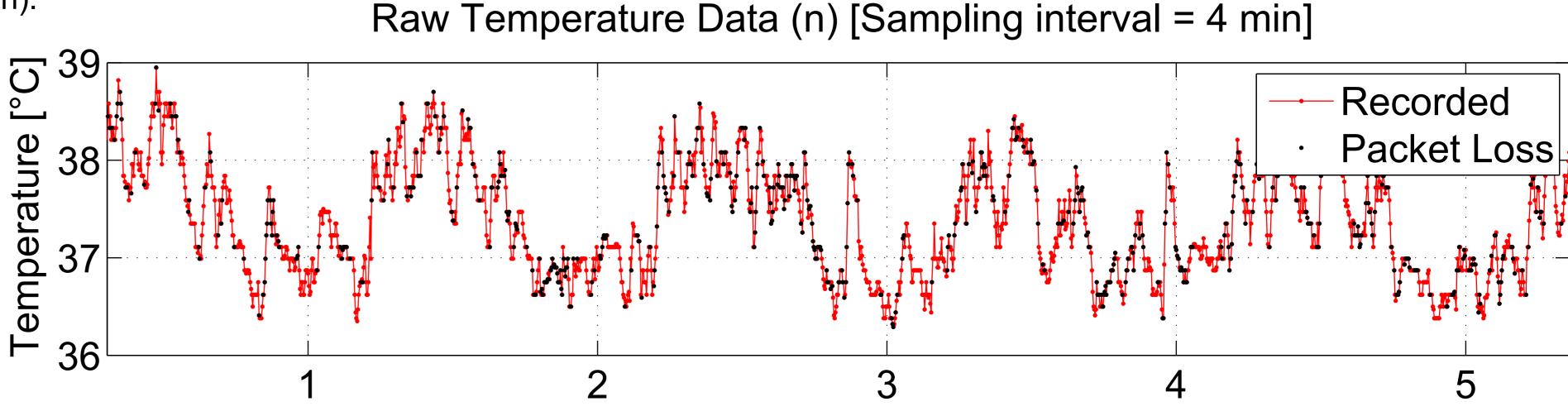



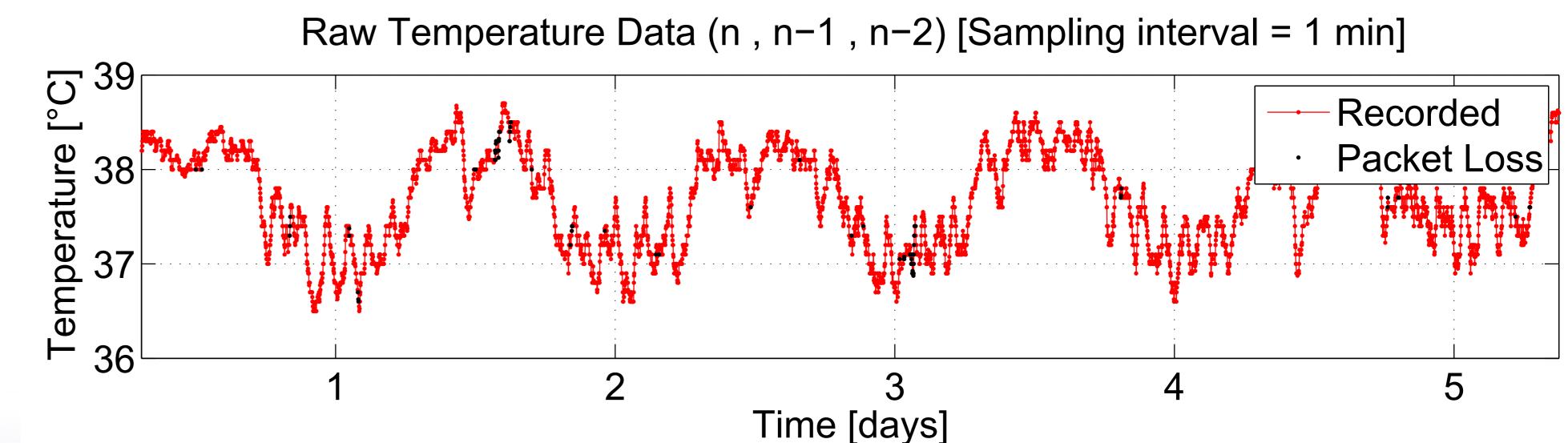

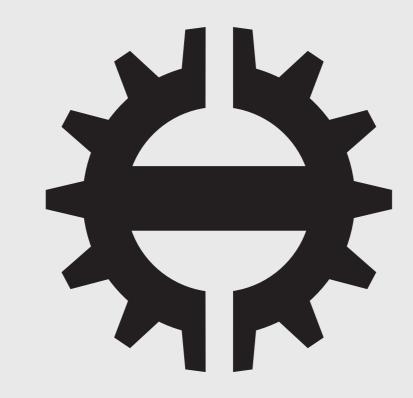

TAMPERE UNIVERSITY OF TECHNOLOGY

